

Instructions of Installing Temperature Controller

1. How to replace temperature controller:

a. At the bottom of the temperature controller there is a lock clip for locking on the temperature controller (see Fig.1).

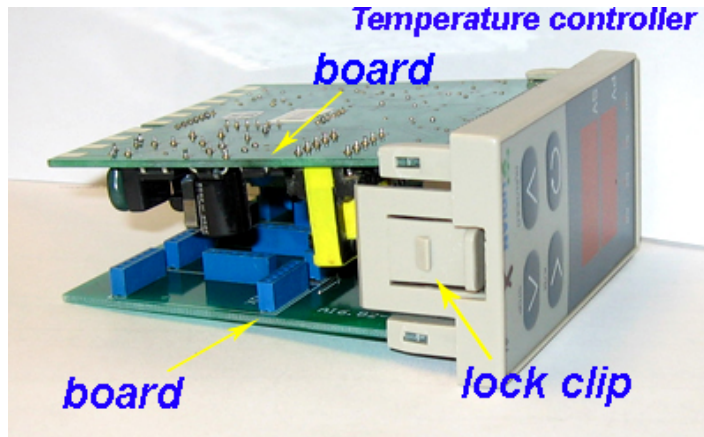


Fig. 1 the view of temperature controller

b. For taking the temperature controller out from the control panel at the bottom of the furnace, please push the lock clip at the bottom of the temperature controller by a screw driver to release the lock (see Fig.2).



Fig.2 uninstallation of temperature controller, step 1

c. Gently pull out the temperature controller by hand while the lock released, shown as Fig.3.

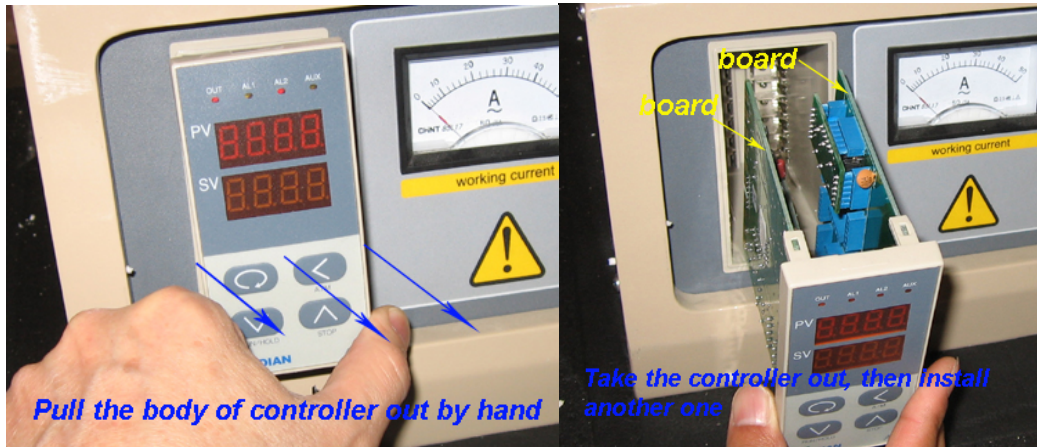


Fig.3 uninstallation of temperature controller, step 2 and step 3

d. Install a new temperature controller by putting it into the panel and slide in directly.

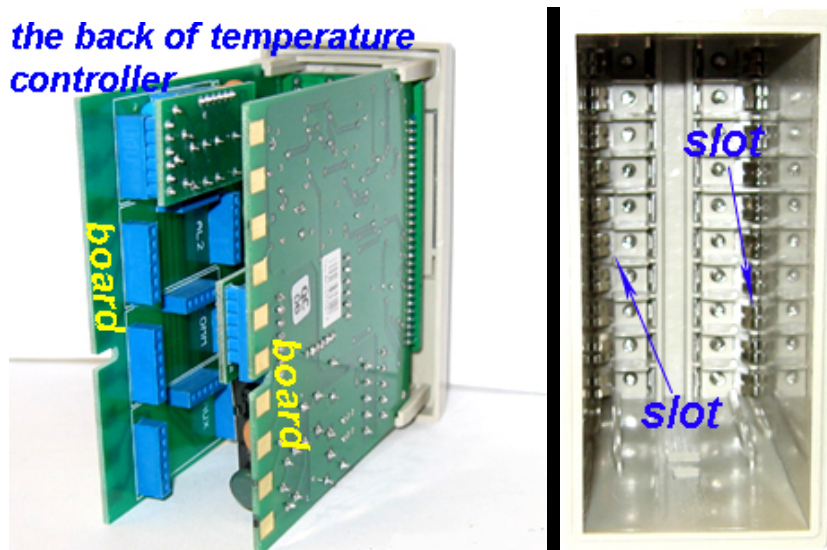


Fig. 4 the view of back of temperature controller and slots

e. Please make sure if two ends of the boards insert into two slots exactly and push the temperature controller all way down to lock it into the panel (Fig. 4). Then run your furnace again.



MTI Corporation

2700 Rydin Road, Unit D, Richmond, CA 94804, USA

Tel: 510-525-3070

Fax: 510-525-4705

E-mail: info@mtixtl.com

Website: www.mtixtl.com