

Trouble shooting

GENERAL OPERATIONS MANUAL



860 South 19th Street, Richmond, CA 94804, USA
TEL: (510) 525 – 3070 FAX: (510) 525 – 4705
Website: www.mtixtl.com Email: info@mtixtl.com

Contents


Troubleshooting 1, Vacuum Problem.....2

Troubleshooting 1, Vacuum Problem

❖ For not able to vacuum the chamber,

First, loosen the Hex Nut.

Second, secure the lock set(towards inside). This will help secure the chamber door.

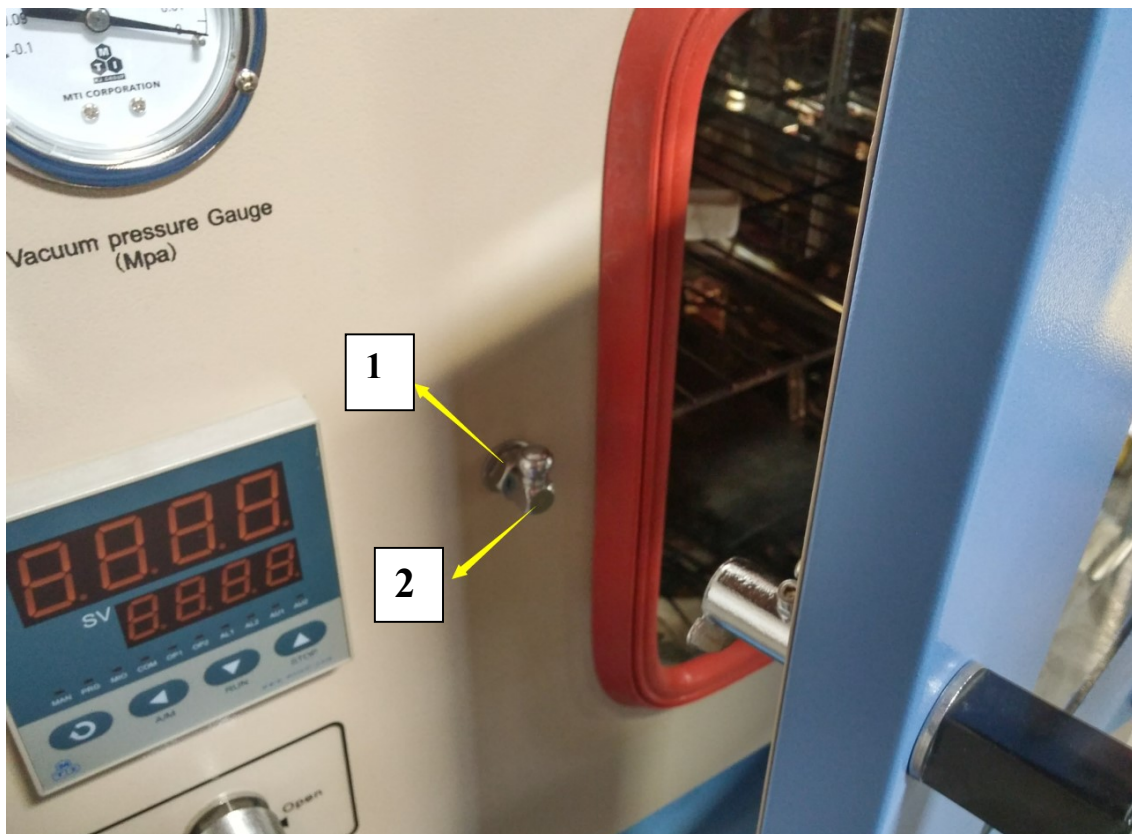

❖ If it did not working, please add some nut gasket at the glass rack (located inside the door.)

