


Coolant Circulating Tank


Structure

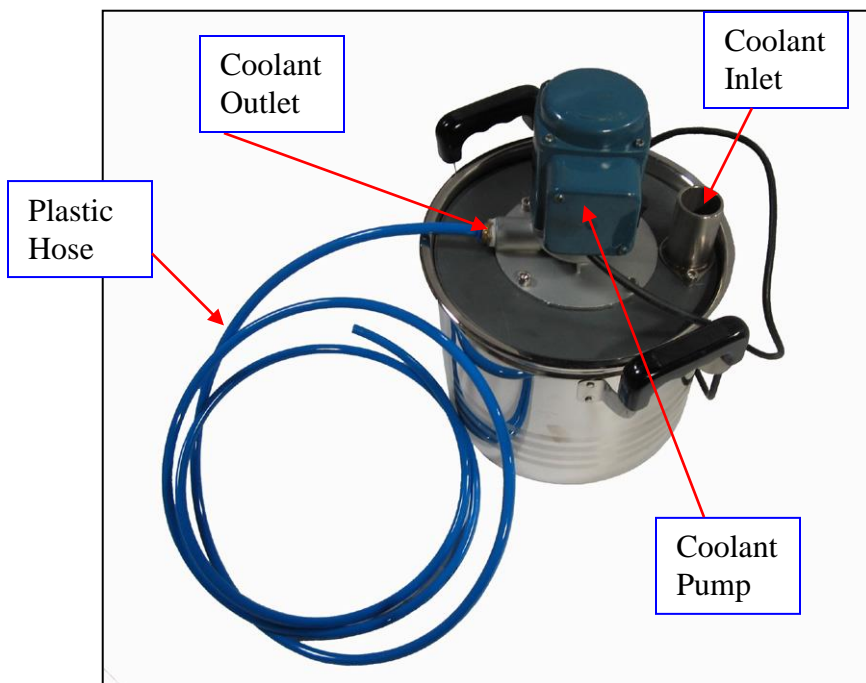

Operation

1. Mix the coolant with water according to the required mixing proportion, and then pour it to the tank.
2. Connect **coolant inlet** to the **drain of device** needs to be cooled down, with a **drain hose** having Outer Diameter of 1". (MTI will not provide this drain hose because it can be easily found in store)
3. Mount one **end of plastic hose** to the **coolant outlet**, and **another end** to the device needs to be cooled down.
4. Power on the machine to start cooling.
5. When finished, press the button to uninstall the plastic hose from the coolant outlet.

Push this button to release the plastic hose from coolant inlet

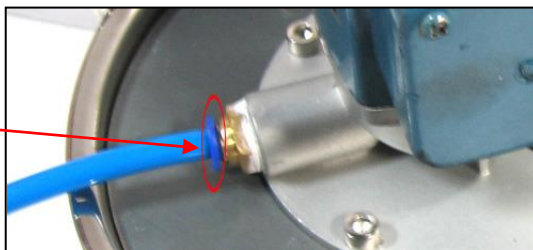

860 South 19th Street, Richmond, CA 94804, USA

Tel: 1-510-525-3070 Fax: 1-510-525-4705

Web: www.mtixtl.com Email: info@mtixtl.com